



Product Manual

SAVIEW[®] (mouse/rabbit-AP) IHC kit

Catalog #: ADI-950-120

(For Mouse and Rabbit Primary Antibodies)



Product Manual

USE FOR RESEARCH PURPOSES ONLY

Unless otherwise specified expressly on the packaging, all products sold hereunder are intended for and may be used for research purposes only and may not be used for food, drug, cosmetic or household use or for the diagnosis or treatment of human beings. Purchase does not include any right or license to use, develop or otherwise exploit these products commercially. Any commercial use, development or exploitation of these products or development using these products without the express written authorization of Enzo Life Sciences, Inc. is strictly prohibited. Buyer assumes all risk and liability for the use and/or results obtained by the use of the products covered by this invoice whether used singularly or in combination with other products.

LIMITED WARRANTY; DISCLAIMER OF WARRANTIES

These products are offered under a limited warranty. The products are guaranteed to meet all appropriate specifications described in the package insert at the time of shipment. Enzo Life Sciences' sole obligation is to replace the product to the extent of the purchasing price. All claims must be made to Enzo Life Sciences, Inc., within five (5) days of receipt of order. THIS WARRANTY IS EXPRESSLY IN LIEU OF ANY OTHER WARRANTIES OR LIABILITIES, EXPRESS OR IMPLIED, INCLUDING WARRANTIES OF MERCHANTABILITY, FITNESS FOR A PARTICULAR PURPOSE, AND NON- INFRINGEMENT OF THE PATENT OR OTHER INTELLECTUAL PROPERTY RIGHTS OF OTHERS, AND ALL SUCH WARRANTIES (AND ANY OTHER WARRANTIES IMPLIED BY LAW) ARE EXPRESSLY DISCLAIMED.

TRADEMARKS AND PATENTS

Several Enzo Life Sciences products and product applications are covered by US and foreign patents and patents pending. Enzo is a trademark of Enzo Life Sciences, Inc.

**FOR RESEARCH USE ONLY.
NOT FOR USE IN DIAGNOSTIC PROCEDURES.**

TABLE OF CONTENTS



Please read entire booklet before proceeding with the assay.



Carefully note the handling and storage conditions of each kit component.



Please contact Enzo Life Sciences Technical Support if necessary.

Intended Use.....	2
Test Principle	2
Introduction	3
Reagents Supplied.....	3
Precautions	4
Storage	4
Other Materials Needed	4
Specimen Preparation.....	4
Reagent Preparation	5
References.....	8
Contact Information.....	10

INTENDED USE

This product is intended for qualitative immunohistochemistry with normal and neoplastic formalin-fixed, paraffin-embedded tissue sections, to be viewed by light microscopy.

This kit is designed to label specific primary antibodies immunohistochemically on tissue sections. The immunohistochemical protocol defined in this brochure is just a guideline. We encourage the individual laboratory to optimize its own protocol based upon tissue fixation conditions, primary antibodies employed, and the user's experience. These reagents were tested and quality controlled using tissue sections, however, they can also be optimized for cell smears and cytospin preparations. Sufficient reagents are provided to run 50 to 100 tests.

TEST PRINCIPLE

The high affinity of the non-covalent interaction between biotin and streptavidin (1×10^{15}) forms the basis for this immunostaining kit. It requires the formation of an irreversible and specific linkage between biological macromolecules. The immunohistochemical applications of the interaction between avidin and biotin were introduced by Bayer et al. (1979), who described techniques for generating active biotinyl compounds such as biotin-N-hydroxysuccinimide and biotin hydrazine and for conjugating them to various organic compounds, including immunoglobulins and alkaline phosphatase (ALP). Streptavidin (SA) is a tetrameric protein (mol. wt. $4 \times 15,000$), isolated from the *actinobacterium Streptomyces avidinii* (Chaiet & wolf, 1964). Streptavidin can bind to four molecules of biotin. Streptavidin gives superior results compared to avidin because its isoelectric point is closer to neutral pH, whereas avidin is positively charged at physiological pH. Streptavidin does not carry any carbohydrate side chain, whereas avidin is composed of 70% of carbohydrate. Because of this, SA does not have the tendency to bind non-specifically. Primary antibodies bind to the target antigens in the tissue sections and conjugated secondary antibodies bind specifically to these receptor antibodies. A biotin-conjugated secondary antibody, in turn, is traced by a streptavidin-conjugated enzyme and can be visualized by an appropriate substrate/chromogen.

INTRODUCTION

Immunohistochemical techniques are spreading rapidly and the practice of anatomic pathology has undergone a revolutionary change since the development of these procedures (Nadji & Morales, 1983). Because of their versatility, sensitivity, and specificity, immunoperoxidase stains are invariably the best stains when appropriate antibodies are available. With the ever increasing number of antibodies against cellular antigens, immunoperoxidase techniques now provide a powerful tool to resolve a wide array of diagnostic pathology. All immunohistochemical techniques require the specific antibody employed to be labeled that they can be easily seen when attached to cellular antigens. At the same time the sensitivity of the immunoperoxidase techniques are central to wide variety of specific antigen localization. Alkaline phosphatase based kits are of special value for staining tissues that have high endogenous peroxidase activity. Alkaline phosphatase is unaffected by endogenous peroxidase and therefore results in cleaner background. Our kit is based on direct SA-AP conjugate technology. The linker reagent is a cocktail of biotinylated anti-mouse and anti-rabbit immunoglobulins, capable of labeling mouse and rabbit primary antibodies.

REAGENTS SUPPLIED

- Bottle 1 Alkaline Phosphatase Enhancer: 10 mL clear solution. Used to increase the intensity of signal.
- Bottle 2 Linker Reagent: 10 mL clear yellow solution of biotinylated anti-mouse and anti-rabbit immunoglobulins.
- Bottle 3 ALP Tracer Reagent: 10 mL clear cherry red solution of conjugated streptavidin alkaline phosphatase.

Detection Reagents:

- HIGHDEF® red IHC chromogen (AP) -- Buffer: 15 mL
- HIGHDEF® red IHC chromogen (AP) -- Chromogen: 1 mL
- One empty mixing bottle

PRECAUTIONS

- i) Several reagents in this kit contain sodium azide. Follow instruction provided by your local authorities for disposal. If disposed in the sink, flush the drain pipe to avoid a reaction of sodium azide in the plumbing system.
- ii) Once you start the immunostaining process, do not let tissue sections dry because it can cause undesirable background staining and artifacts.
- iii) Interpretation of the results will be the sole responsibility of the user.

STORAGE

All of the kit components should be stored at 2-8°C. Do not freeze. Do not use beyond the expiration date stated on the label.

OTHER MATERIALS NEEDED

- Xylene or dewaxing agents
- Absolute alcohol
- Distilled or deionized water
- Counterstain
- IHC wash buffer (Prod. no. ADI-950-235)
- HIGHDEF[®] IHC mount (Prod. no. ADI-950-261)
- Oven or incubator
- Staining jars
- Microscope slide
- Microscope
- Diamond pencil
- Absorbent pads
- Primary antibodies

SPECIMEN PREPARATION

The purpose of immunohistochemistry is to localize, identify, and (whenever possible) quantitate cell and tissue constituents. Optimal fixation of the tissue is the key to good immunostaining. Successful immunostaining depends on the accurate preservation of the epitopes looked for in its structural context. We recommend our users to read an excellent article on tissue preparation by Larsson (1993) to understand the need and importance of optimal tissue fixation.

REAGENT PREPARATION

Except for the chromogen, all reagents are provided in ready-to-use form.

Preparation of Substrate/Chromogen

Aliquot 1 mL of HIGHDEF® red IHC chromogen (AP) buffer in a mixing bottle. Add one drop (~20 µL) of HIGHDEF® red IHC chromogen (AP) chromogen. Replace tip, mix, and allow solution to reach room temperature before using.

Note: The working chromogen-substrate solution should be prepared fresh and used within 20-30 minutes of preparation. Any solution not used during this period should be discarded.

Positive and Negative controls

Each immunostaining run should include a known positive and a negative control to assure the proper functioning of staining system and valid interpretation of the results.

Positive control:

A tissue which is known to contain the desired antigen and has given positive staining in the past.

Negative control: One of the following should be used as negative control:

- i) Instead of primary antibody, use the normal non-immune serum from the same species of animal in which the primary antibody was raised.
- ii) Use buffer in which the primary antibody was diluted.
- iii) Use a tissue known not to contain the desired antigen.
- iv) Absorb the primary antibody with the appropriate antigen and use it instead of primary antibody.

Staining Protocol

- Step I Removal of paraffin wax: Deparaffinize tissue sections according to the established procedure in your lab and bring tissues to wash buffer.
- Step II Endogenous Alkaline Phosphatase Blocking: This is an optional step and should be performed only if tissue is suspected to have high endogenous alkaline phosphatase activity. Apply enough drops of Alkaline Phosphatase Blocker to cover tissue. Incubate for 5-10 minutes at room temperature.
- Step III Washing: Drain off excess reagent. Rinse with wash buffer three times for 1 minute each time. Drain off excess buffer and carefully wipe slide around the tissue to remove excess buffer from the glass leaving the tissue wet.

Immunostaining Protocol I

This protocol is recommended for optimally fixed tissues with abundant antigens in the tissue and for high affinity primary antibodies.

- Step I Primary antibody: Apply enough drops of primary antibody to cover the tissue section. Incubate according to the manufacture's recommended conditions. Wash and wipe slides as described above.
- Step II Linker reagent: Apply enough drops of linker reagent to cover the tissue section. Incubate for 10 minutes at room temperature. Wash slides as described previously.
- Step III Tracer Reagent: Apply enough drops of tracer reagent to cover the tissue section. Incubate for 10 minutes at room temperature. Wash slides as described previously.
- Step IV Substrate/Chromogen: Apply working chromogen solution for 5-10 minutes at room temperature for color development. For best results, look under the microscope for signal development. Once desired signal-to-noise ratio is achieved, stop the reaction by washing slides in wash buffer. Note development time and follow it during subsequent incubations.

Immunostaining Protocol II

This protocol is recommended for less than optimally fixed tissues with low antigenic density in the tissue and for low affinity primary antibodies.

- Step I Primary antibody: Apply enough drops of primary antibody to cover the tissue section. Incubate according to the manufacture's recommended conditions. Wash slides as described previously.
- Step II Linker reagent: Apply enough drops of linker reagent to cover the tissue section. Incubate for 20 minutes at room temperature. Wash slides as described previously.
- Step III Tracer Reagent: Apply enough drops of tracer reagent to cover the tissue section. Incubate for 20 minutes at room temperature. Wash slides as described previously.
- Step IV Substrate/Chromogen: Apply working chromogen solution for 5-10 minutes at room temperature for color development. For the best results, look under the microscope for the signal development. Once desired signal-to-noise ratio is achieved, stop the reaction by washing the slide in wash buffer. Note development time and follow it during subsequent incubations.
- Step V Wash slides and counter stain with an appropriate counter stain. Mount and observe staining under the microscope.

Enhanced Staining

If additional signal is desired, treat tissue sections with Alkaline Phosphatase Enhancer for 1 minute after Step III. Drain off excess enhancer and add substrate/chromogen solution without any wash.

REFERENCES

1. Bayer et al., Method Enzymol. 62, 308 (1979).
2. Chaiet & Wolf, Arch. Biochem. Biophys. 106, 1 (1964).
3. Giorno, Diag. Immuno. 2, 161 (1984).
4. Hsu et al., J Histo. Cyto. 29, 577 (1981).
5. Nadji & Morales, Anat. Pathol. 14, 767 (1983).
6. Nagle et al. J Histo. Chem. 31, 1010 (1983).
7. Larsson, Applied Immunohistochem. 1, 2, (1993).
8. Petrusz & Ordronneau, In Polak & Van noorden eds., Immunocytochemistry: Practical applications in pathology and biology. Bristol Wright-PSG, 212, (1983).



Product Manual

NOTES



Product Manual

www.enzolifesciences.com

Enabling Discovery in Life Science®

GLOBAL HEADQUARTERS

Enzo Life Sciences Inc.
10 Executive Boulevard
Farmingdale, NY 11735
Toll-Free: 1.800.942.0430
Phone: 631.694.7070
Fax: 631.694.7501
info-usa@enzolifesciences.com

EUROPE/ASIA

Enzo Life Sciences (ELS) AG
Industriestrasse 17
CH-4415 Lausen
Switzerland
Phone: +41/0 61 926 89 89
Fax: +41/0 61 926 89 79
info-ch@enzolifesciences.com

For local distributors and detailed product information visit us online:

www.enzolifesciences.com